# User Guide

## NUL-229

English

Español

# NeuLog logger sensors and modules

## NeuLog logger sensors

Welcome to the new age of experimental data acquisition with plug and play smart logger sensors.

NeuLog logger sensors offer a myriad of didactic tools to meet your needs in the field of scientific studies.

The NeuLog system works as a set of independent computer-based modules complete with a data logger, flash memory and a sensor – all in each individual sensor.

Sensors plug into each other and have the ability to connect as a chain of sensors and then stream data to any computer system or smart device.

The NeuLog system can connect to nearly any computer platform or operating system (PC, MAC, iPad, Tablets, Android, smart phones, etc). All that is needed is a USB connection or wireless connectivity.

**Note:** NeuLog products are intended for educational use.

## NeuLog user guide

This is a general guide for all NeuLog sensors and modules instructing how to use them at various operation modes. Specific user guides for each sensor and module can be found at [www.NeuLog.com](http://www.NeuLog.com). Choose "Products" on the main menu, and then choose the relevant sensor or module.

Each specific user guide contains the following sections:

- What is included with the sensor
- Sensor specifications
- Technical background
- Maintenance and storage
- Sensor offset (if needed)

**Note:** There are some sensors (listed below) that due to their specific nature require different instructions than the rest of the NeuLog sensors. For these sensors please make sure to refer to [www.NeuLog.com](http://www.NeuLog.com) before use:

- NUL-209, Photo gate sensor
- NUL-243, GPS

## Quick start procedure

### PC or Mac Computer – Wired Connection

#### Materials needed:

- NeuLog sensor or sensors
- USB-200 USB Module, or WIFI-201/WIFI-202 (using WiFi as USB connection)
- USB to mini USB cable (included with the USB-200 module)


Your sensor needs to be connected to a USB-200 module. The USB-200 module then connects to a computer via USB (this can also be done with WIFI-201 or WIFI-202). Please note that you cannot plug the sensor directly into the computer.

The browser-based application can be downloaded for free at [www.NeuLog.com](http://www.NeuLog.com) as well as a full application user guide.

Choose "Downloads" on the main menu and then choose "Software and Application".

## Procedure:

1. Install the NeuLog application.
2. Connect the USB-200 module to the PC or Mac.
3. Connect the sensor or sensors to the USB-200 module (they directly plug together). Please note that no calibration is required for NeuLog sensors. You should see the red LED on the sensor flash as you plug them in.
4. Open the NeuLog application by clicking on the shortcut on the screen.
5. The connected probes are automatically identified by the software; including disconnecting or connecting other probes.

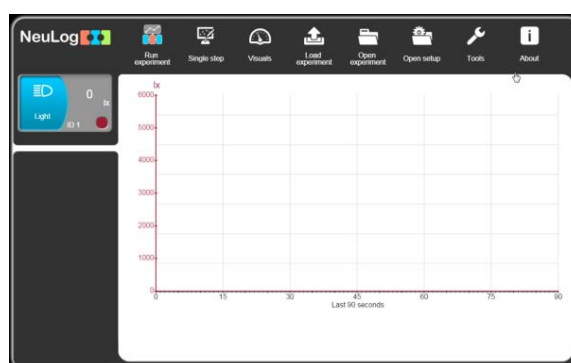6. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

7. Select one of the following function buttons and the software will guide you through the function's options:

- 'Run Experiment' with the connected sensors.
- 'Single Step' experiment.
- 'Visual' display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
- 'Load Experiment' from the sensors' flash memory.
- 'Open Experiment' from a saved file.
- 'Tools' for changing the software settings or sensor ID.

### Tablet, smart phone or any device with WIFI connectivity

#### Materials needed:

- NeuLog sensor or sensors
- WiFi Module (WIFI-201 or WIFI-202)
- Battery Module (BAT-200 or BAT-202)

Your sensor or sensors needs to be connected to a WiFi module. The WiFi module will create a closed NeuLog WiFi network which will stream the NeuLog data to the device of your choosing. Once your device is wirelessly connected to the NeuLog network, you can run experiments and collect data through a browser of your choosing.

**Procedure:**

1. Connect the sensor or sensors directly to the left side of a WiFi module (no wires required).
2. Connect a BAT-200 module to the right side of the WiFi module or BAT-202 to its left side.
3. Although not required, we recommend plugging the BAT-200 to an outlet using a USB to mini USB cord. The WiFi module will run for 60-90 minutes (depending on the sensor) without being plugged in.
4. The WiFi module can be directly powered using a mini USB to USB cord and plugging it into your computer or a wall charger. Please note that this function is only available on WiFi modules with the mini USB port.
5. For further WiFi module instructions or the WIFI-201 or WIFI-202 quick start guide please visit: [www.NeuLog.com](http://www.NeuLog.com). Choose "Products" on the main menu, then choose "WiFi Communication module".
6. The WiFi module indicator lights will flash. Take no action until the LED turns blue. This can take up to one minute.
7. Take your tablet or smart phone, go to the WiFi settings and select the NeuLog network (NeuLog0184 for example) which matches the ID found on the back of the WiFi module (0184 for this example).
8. Give your device about 20 seconds to connect to the WiFi module.

9. Once the device is connected, go to your browser and type the website **wifi201.com** into the URL bar, then wait for 30-60 seconds.
10. The connected probes are automatically identified by the software including disconnecting or connecting other probes.
11. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.
12. If you are not the first one to connect to this WiFi unit, enter viewer mode and view the screens of the user master who was the first to connect.
13. If you are the first one to connect (the Master), you can select one of the following function buttons and the program will guide you through the function's options:
  - a. **'Run Experiment'** with the connected sensors.
  - b. **'Single Step'** experiment.
  - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
  - d. **'Load Experiment'** from the sensors' flash memory.
  - e. **'Open Experiment'** from a saved file.
  - f. **'Tools'** for changing the program settings or sensor ID.

14. You can click on the 'Release Master' button in the 'Tools' options and all the connected viewers will have a new 'Be Master' button. The first one to click becomes the 'Master'.
15. The 'Tools' also includes a 'WiFi Options' button that enables to connect the WiFi module as a client to the WiFi network and to connect to the module through this network. This option allows using the internet while using the NeuLog sensors.
16. If you have a newer generation of WIFI-201 module (with the USB icon on the cover) or a WIFI-202 module this can be used as a USB-200, meaning you can tether the probes directly to the computer using the WiFi module. To engage operation in this fashion, press three times on the panel button. Repeating this operation will return the module to WiFi mode.

### Tablet, smart phone with Bluetooth connectivity

#### Materials needed:

- NeuLog sensor or sensors
- BLT-200 – Bluetooth Module (requires IOS 10 and up for Apple devices and BLE (Bluetooth Low Energy) for Androids devices)

Download the NeuLog app from the "App Store" or from "Google Play".

**Procedure:**

1. Connect the sensor or sensors directly to the right side of the Bluetooth module (no wires required).
2. Turn the Bluetooth module ON by pressing the BLT-200 button, the blue LED will flash.
3. Run the app. The app will search for the module and connect automatically. After connecting, the blue LED on the module will light constantly, and at the app the Bluetooth sign (on the right top corner) will turn green.
4. If the connection fails, press the gray Bluetooth sign (on the right top corner) to refresh the connectivity.
5. If there is more than one Bluetooth module available, the app will ask you to choose which module to connect.
6. The connected sensors are automatically identified by the software including disconnecting or connecting other probes.
7. Once a sensor module box appears on the left side of the screen, the sensor will give a live reading in its module box and on the graph appearing on the screen and you can begin experimentation.

8. You can select one of the following function buttons and the software will guide you through the function's options:
- a. **'Run Experiment'** with the connected sensors.
  - b. **'Single Step'** experiment.
  - c. **'Visual'** display with four options for one of the connected sensors values: Digital, Bar, Analog, gauge and Accumulated graph.
  - d. **'Load Experiment'** from the sensors' flash memory.
  - e. **'Open Experiment'** from a saved file.
  - f. **'Tools'** for changing the program settings or sensor ID.

### Operation with NeuLog Graphic Display Module (viewer)

#### Materials needed:

- NeuLog sensor or sensors
- VIEW-101 Graphic Display Module
- BAT-200 or BAT-202 Battery Module

**Procedure:**

1. Connect the sensor or sensors to the viewer.
2. Connect the BAT-200 or the BAT-202 to the viewer.
3. The VIEW-101 will automatically detect the sensor and you will see it appear on the left side of the screen.
4. Once the sensor appears it will be monitoring data in real time.
5. To run an experiment and collect data, press the run icon (little green person).

**Off-line experiments**

Off-line experiments are when you do not have a sensor connected directly to a computer, tablet, smart phone, or NeuLog viewer.

**Materials needed:**

- NeuLog sensor or sensors
- BAT-200 or BAT-202 Battery Module

**Materials needed to configure your offline experiment:**

- USB-200 Module, WiFi module (WIFI-201 or WIFI-202) or VIEW-101 Module
- USB to mini USB cable (included with the USB-200 module).

**Procedure:**

1. To determine the experiment settings, first connect to your chosen device (PC, Mac, tablet, smart device) and wait until the sensor search has ended.
2. Click on each sensor's module box, on the left side of the screen, and set the experiment parameters (Range, Duration, Rate and Trigger) for this sensor. The setting parameters will be saved in the sensor's flash memory.
3. Connect the sensor or sensors to a charged BAT-200 or BAT-202 module.
4. When ready to collect data, press the "Start/Stop" button on the sensor. A red light will turn on for the duration of the experiment.
5. After the experiment has concluded (the "Start/Stop" button is pressed again or the data collection period ends). Connect the sensor or sensors to your chosen device (you can find full instructions in each device's quick start procedure section).
6. Open the NeuLog application.
7. Click the 'Load experiment' icon.
8. Select which experimental data to upload (up to 5 experimental runs can be stored at one time on the sensor).

## Operation with NeuLog Digital Display Module (viewer)

### Materials needed:

- NeuLog sensor or sensors
- VIEW-200 Digital Display Module
- BAT-200 or BAT-202 Battery Module

Prior to data collection, it is recommended that you configure your sensors' settings to your specific experimental needs. Instructions for off-line experiments configuration procedure can be found in the "Off-line experiments" section in the previous page.

If you only use the VIEW-200 to take live readings without collecting data it is not required to configure your sensors' settings (unless, if needed, for changing the sensors' range).

**Procedure:**

1. Connect up to 5 NeuLog sensors of your choice together and into the left side of your viewer.
2. Connect a charged battery module into the right side of the viewer. The screen will automatically turn on.
3. The viewer will begin searching for sensors at this time – the screen will read "Search... %". This process should take roughly 25 seconds to complete.

**Note:** If no sensors are found, disconnect and reconnect the battery module.

4. Once the sensors have been detected, you can begin collecting data directly onto the internal memory of each sensor. To do this you simply need to press the "Start/Stop" button on the front of each sensor you wish to collect data with.
5. You can change which sensor data is displayed on the VIEW-200 by pressing the button on the front of the module.

# NeuLog Soil Moisture Logger Sensor Guide NUL-229

## NeuLog soil moisture logger sensor NUL-229

The NeuLog soil moisture sensor can be used for any science experiment which utilizes accurate soil moisture levels such as in the fields of Environmental Science, Ecology, Animal Behavior, Biology, Chemistry, etc.

The sensor comes pre-calibrated so you can start experimentation right out of the box using this guide.

Among hundreds of possible experimental subjects that can be studied with the NUL-229 sensor are: monitoring habitats, ecosystem health, animal behavior, soil chemistry, soil comparisons, ecological monitoring, and many more.

This logger sensor is based on measurements of the vacuum pressure in a tensiometer. A tensiometer is a closed tube full with water with a special ceramic part in its end. In dry soil, water goes out by diffusion through the ceramic holes and vacuum pressure is created in the tensiometer. The vacuum pressure is proportional to the soil dryness.

The soil moisture sensor's measurement units are:

- Centibar (cbar): The bar is a non-SI unit of pressure defined as 100,000 Pascals. The centibar is defined as one kilopascal.
- Kilopascal (kPa): the SI derived unit of pressure.
- Percentage (%)

## Setting up the soil moisture probe

The NeuLog soil moisture sensor requires only a minimal setup prior to each use. Be sure to properly configure your NeuLog sensor using one of the following guides (computer/smart device/viewer) prior to data collection.

1. Remove the soil moisture probe's cap which is wired directly into the sensor's body.
2. Immerse the ceramic cone in tap water. It is recommended to leave it in the water for 15 minutes.
3. Fill the probe's ceramic cone with tap water.

**Note:** More water will last longer and provide more accurate results though be sure to allow some room for air at the top of the column.

4. Replace the plastic cap and firmly press down to form a strong seal.
5. Gently press the probe (ceramic cone side first) vertically into the soil that you want to test.
6. Ensure that the entire ceramic cone has been pressed into the soil.
7. Begin collecting data.

### Included with sensor

- NeuLog General Guide.
- Blumat soil moisture probe attached directly to the NeuLog sensor's body.

Sensors specifications			
	cbar	kPa	% moisture
Range and operation modes	-20 to 50 cbar	-20 to 50 kPa	0-100%
ADC resolution	15 bit		
Resolution	0.01 cbar	0.01 kPa	5%
Max sample rate (S/sec)	100		

**Experiment Duration:** 1 second to 31 days.

### Sensor's features

- Fully digital data.
- Rugged plastic ergonomic case.
- Pushbutton switch for Start/Stop experiments in off line mode.
- LED indicator of experiment status (blinks while collecting data).
- Pre-calibrated sensing equipment.

**Note:** NeuLog products are intended for educational use.

## Videos and experiment examples

- Videos, literature and other probes can be found at [www.neulog.com](http://www.neulog.com).
- In order to access the soil moisture sensor's page, choose "Products" on the main menu and then "Soil Moisture Logger Sensor".
- In order to access the soil moisture sensor's experiments, choose "Example Labs":

Soil Moisture (B-20)

## Technical background

The philosophy behind NeuLog's plug and play technology is based on each sensor's ability to store its own data due to an internal flash memory chip and micro-controller in each plastic NeuLog body. This technology allows the sensor to collect and then store the digital data in the correct scientific units (for example: °C, °F, Lux, %, ppm).

The sensor is pre-calibrated at the factory. The built-in software in the logger can be upgraded for free at any time using the provided firmware update.

The soil moisture sensor is a very technological complex system which utilizes a Blumat soil moisture probe in conjunction with the NeuLog sensor's hardware and software.

After the probe has been filled with water and placed into a soil sample, the specialized ceramic cone tip allows for water diffusion across a gradient. A vacuum is formed inside the probe when the cap is sealed and the sensor is automatically calibrated.

When the soil has less moisture than the probe, water diffuses out of the ceramic cone creating a larger volume vacuum. When the soil's moisture content is higher than the probe's, water will diffuse into the probe through the ceramic cone creating a smaller volume vacuum.

The NeuLog soil moisture sensor directly measures the pressure of the vacuum so the soil moisture reading is presented in pressure units.

## Maintenance and storage

- Never submerge the NeuLog plastic body in any liquid.
- Do not allow liquid into the soil moisture sensor's body.
- After use, gently wipe away any foreign material from the soil moisture sensor.
- Store in a box at room temperature out of direct sunlight.

## Warranty

We promise to deliver our module free of defects in materials and workmanship for a period of 3 years from the date of purchase. Our warranty does not cover damage of the product caused by improper use, abuse, or incorrect storage. Sensors with a shelf life such as ion selective probes have a warranty of 1 year. Should you need to act upon the warranty please contact your distributor. Your module will be repaired or replaced.

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